

## Immunofluorescence- **KRT14**, **MUC5B**, and **MUC5AC**

**Purpose:** Stain slides of 5µm paraffin sections from the HTC of adult human lung for **KRT14**, **MUC5B**, and **MUC5AC** with antigen retrieval.

### Day 1

1. Paraffin sections are placed at 60°C for 2 hours to overnight to melt paraffin.
2. Paraffin sections are then placed in xylene 3X for 10 minutes each, followed by 3X in 100% ETOH for 3 minutes each, 95% ETOH for 3 minutes each, and 70% ETOH for 3 minutes each. Slides are then placed in 1X PBS for 5 minutes to completely rehydrate tissue.
3. Antigen retrieval, pH 6.0 (times will vary according to microwave).
4. 10 mM sodium citrate, pH 6.0, and heat in a microwave at 96°C. \*We usually do a series of three runs (6-7 minutes each run) to equal the time/temp because of evaporation (refill coplin jars with dH<sub>2</sub>O).
5. Microwave according to instructions on microwave.
6. Cool on countertop, 15 min.
7. Rinse with dH<sub>2</sub>O
8. 1X PBS, 5 min.
9. Block in 4% Goat serum/PBS-T, 2 hours at RT.
10. For mouse anti-**KRT14** (**AB7800**, **Abcam**) dilute 1:100, for rabbit anti-**MUC5B** (**SC-20119**, Lot # B1003 Santa Cruz) Dilute 1:100, for mouse anti-**MUC5AC** (Abcam, AB3649, Lot# GR135303-1) dilute 1:200 in blocking buffer. Spin down in µfuge for 10 minutes and incubate on tissue overnight @ 4°C.

### Day 2

1. Rinse slides in PBS-T 3X, 5 min.
2. Apply secondary antibody, Goat Alexa Fluor 488 anti-mouse IgG3 (**A21151**, Lot# 1563682, for anti-**KRT14**) at 1:200, Goat Alexa Fluor 555 anti-rabbit IgG (**A21428**, Lot# 1608466 for anti-**MUC5B**) at 1:200, Goat Alexa Fluor 647 anti-mouse IgG1 (**A21240**, Lot# 1562685, anti-**MUC5AC**) at 1:200, in blocking buffer. Spin down in µfuge for 10 min, apply to tissue and incubated at room temperature for 1 hour.
3. Rinse in PBS-T 3X, 5 min.
4. Dilute DAPI 1:2000 and apply to slides for 10 min.
5. Wash in PBS-T 3X, 5 min.
6. Rinse slides in 0.1M PB or TB, 3X, 5 min.
7. Add 1 drop of Prolong Gold anti-fade mounting medium (P36930).
8. Coverslip with Gold Seal Coverslip (Cat# 3422 Electron Microscopy Sciences, 22 X 22 mm).
9. Allow Prolong Gold to cure overnight at room temp in light sealed box.

**Tissue Used:**

**LMH-15-D007-RLL-10B2.17**

**Gender: Male**

**Age: 4 Years**

## **Appendix**

### **Antigen Retrieval Solution**

18ml Solution A- 0.1M Citric Acid (Sigma, C1909) pH 2.5

82ml Reagent B- 0.1M Sodium Citrate (Sigma, S4641) pH 8.2

1L dH<sub>2</sub>O

pH 6.0